


ProPac 3R SCX , MAbPac SCX, ProPac Elite WCX 色谱柱上盐梯度下分析三抗

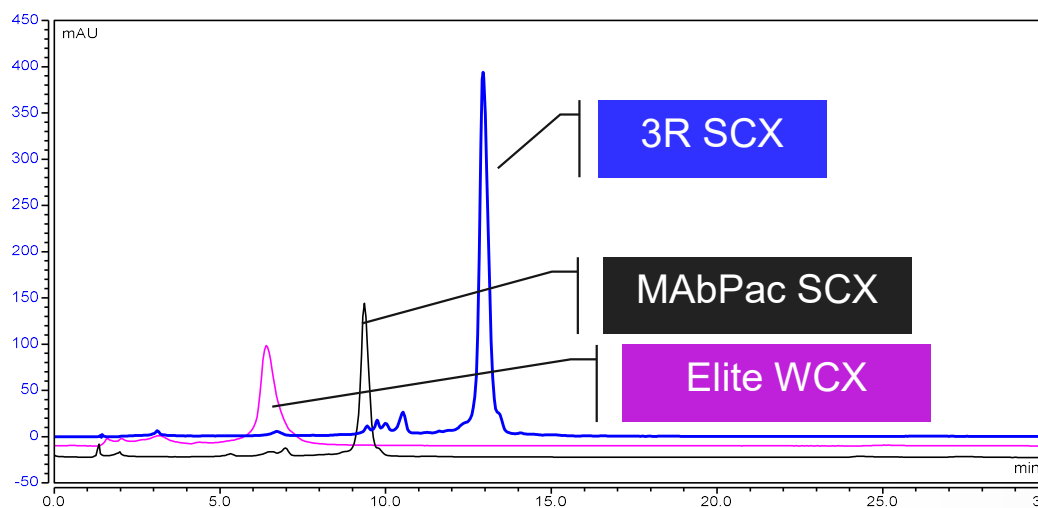

Column: ProPac 3R SCX, 3 μ m
Format: 4x100 mm
Buffer A: 20mmol MES pH 6.7
Buffer B: 20mmol MES +500 mmol NaCl pH 6.7
Flow rate: 0.32mL/min/0.2mL/min
Gradient: 14min(25%-60%)/44min(25%-60%)
Temperature: 30°C
Detection: 280nm

- 三抗电荷异构体分析, 对比0.3ml/min和0.2ml/min的流速下, 20分钟梯度和50分钟梯度, 50分钟可以将难分离的碱性峰分得更好, 流速对分离结果影响不大, 降低流速没有很严重的扩散效应-----省溶剂

